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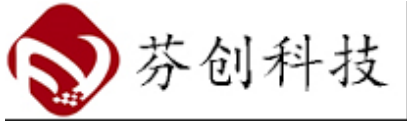
Observation unit with polarized illumination

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I. Overview

1. Features

- 1). Polarized field observation is possible with polarized illumination. The
brightness of the illumination light can be adjusted by the aperture
stop.
- 2). Compatible with infinity correction objective lenses for the visible
range. The objective lens mounting screw is φ 26 x 0.706.
- 3). Built-in tube lens with a focal length of 200 mm.
- 4). C-mount camera, halogen lighting, and LED lighting can be installed.
- 5). Various filter can be attached.
- Note): Objective lens, C-mount camera, lighting, and filter are sold
separately.

2. Specification

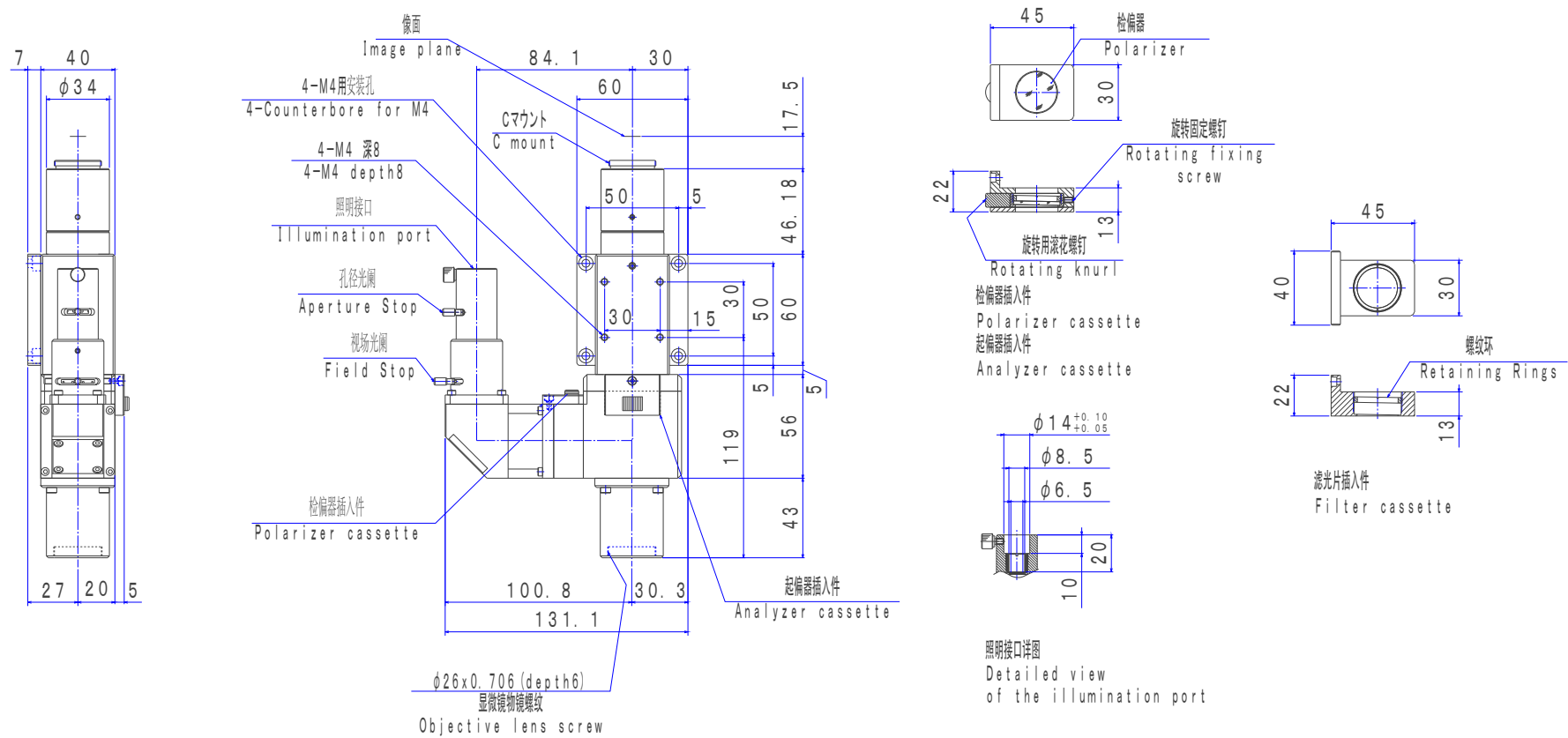
- Product number: OUCI-3-PFU
- Tube lens focal length: 200mm
- Observation image: bright field (Polarized observation)
- Camera mounting direction: vertical
- Camera mount: C mount
- Compatible camera size: 1.2 type (diagonal 21.4 mm) or less
- Lighting: Coaxial epi-illumination
- Lighting mounting shape: Please refer to the outline drawing.
- Corresponding objective lens screw: φ 26 x 0.706
- Corresponding objective lens: Infinity correction objective lens for visible
range
- Mass: 950g
- Polarizing filter : Polarizer has a rotation mechanism.
Analyzer has a rotation mechanism.
- Filter : φ 25.4mm、25.0mm t≤8.5mm (With 2 filter cassettes)

II. Precautions

- 1). Do not use it for any purpose other than its intended use as a
microscope.
- 2). Please note that the magnification will be different if you use an
objective lens designed with a tube lens other than f = 200mm.
- 3). A finite objective lens cannot be used.

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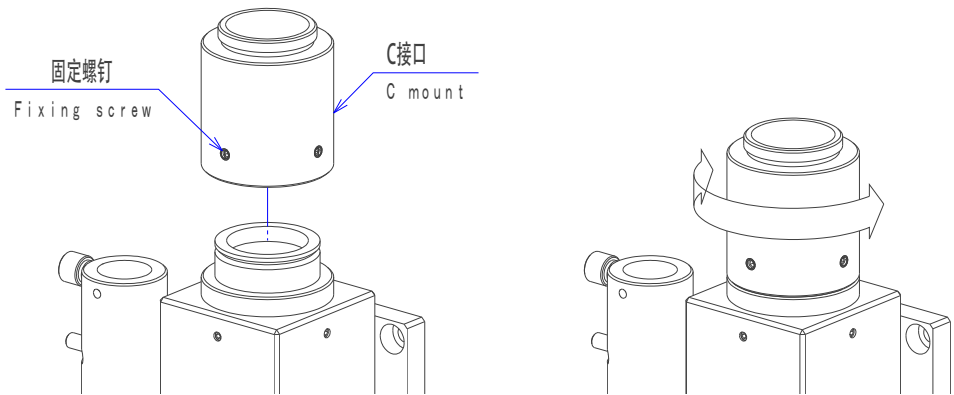
IV. Handling method

1. Installation of OUCI-3-PFU

Fix OUCI-3-PFU with M4 screws. (There is a counterbore hole for the M4 hexagon socket head cap screw.)


2. Installation of camera (sold separately)

Loosen the fixing screws (2 places) and remove the C mount. After attaching the camera to the C mount, put it back. Adjust the rotation of the camera and tighten the fixing screws (2 places).

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
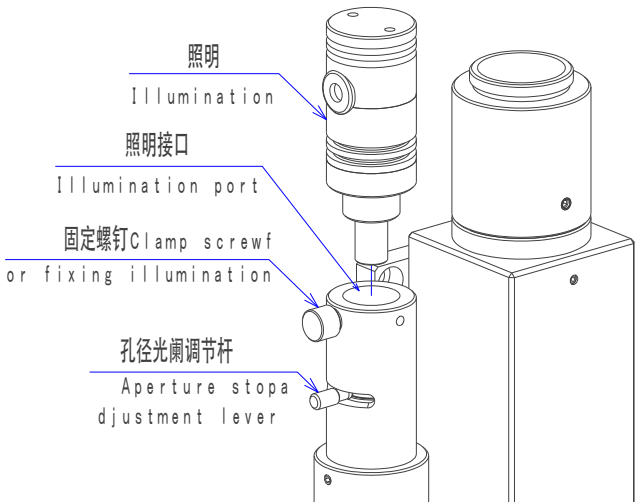
3. Installation of illumination (sold separately)

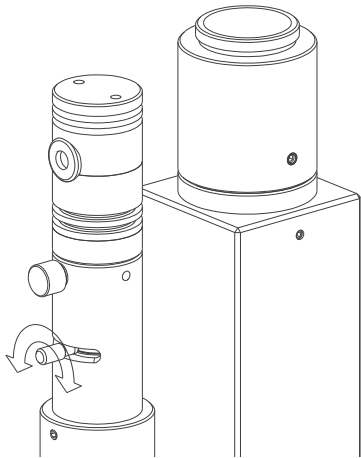Insert the illumination into the illumination port and tighten the fixing clamp screw.
You can adjust the brightness by moving the aperture stop adjustment lever.

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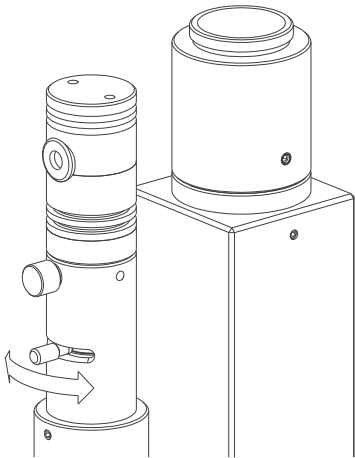Turn clockwise to lock
Turn counterclockwise to release

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4. Installing the filter

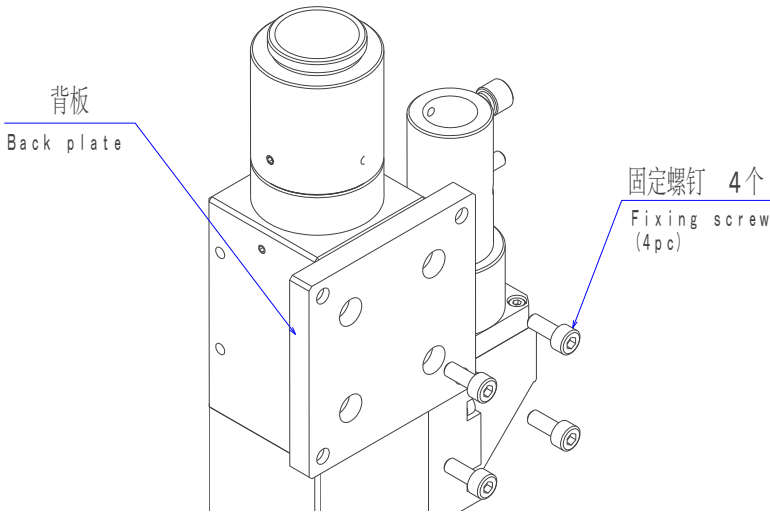
Remove the retaining rings of the attached filter cassette and install the filter.
Remove the polarizer and analyzer cassette attached to the main unit.
Insert the filter cassette in that empty filter slot and tighten it with the fixing screw.
Note) Filter and screw ring spanner are sold separately.

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
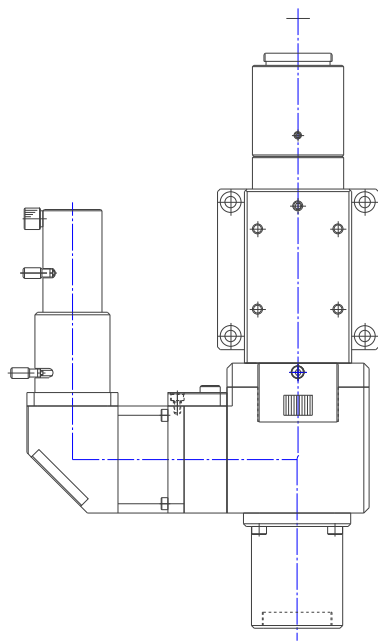
5. Change the mounting orientation

Remove the fixing screw on the back plate. (M4 hexagon socket head cap screw)
Move the removed back plate to the surface you want to change.
Tighten the fixing screw.


出厂时
Factory default

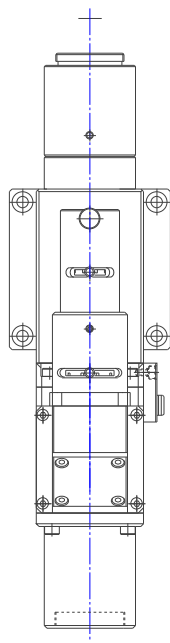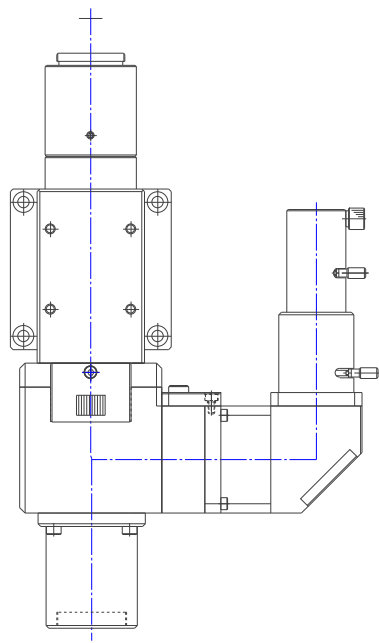安装时 方案1
Change of mounting surface
type 1

安装时 方案2
Change of mounting surface
type 2

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V. Warranty

The warranty period of this product is one year after delivery. In the unlikely event that a malfunction occurs due to our responsibility during the warranty period, please return it to us. We will repair or replace the parts free of charge. However, indirect or consequential damages (including lost profits) caused by the failure of the delivered product are excluded from this warranty. Please refer to our web for details.

VI. Others

- 1). The contents of this manual do not permit the implementation or use of rights such as intellectual property rights owned by SIGMA KOKI Co., Ltd. Reprinting or copying of part or all of this document requires the prior written consent of SIGMA KOKI Co., Ltd.
- 2). The contents of this manual are as of the time of publication and are subject to change without notice.
- 3). The contents of this manual have been carefully prepared, but if you find any mistakes in the description of the materials, please contact the place of purchase or our sales department.
- 4). Combinations other than those described in this manual, or custom specifications may differ from the contents of this manual. If there are products that can be used in combination, also read the instruction manuals for those products.